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لیست محیط کشت های آزمایشگاه میکروبیولوژی مهندسی بهداشت محیط			
Code	Name	Preparation	Abb.
A1	Agar Powder	-	
A2	Azide dextrose broth	Suspend 35g or 70g in 1L of Demin. Water. Dispense into smaller vessels. Autoclave 15 min. at 121C (DO NOT OVERHEAT) Final PH 7.2 ± 0.2 at 25C	
B1	Bile Esculin Agar	Suspend 64.5g in 1L of Demin. Water. Boil to dissolve the medium completely. Autoclave 15 min. at 121C & 15 lbs Final PH 6.6 ± 0.2 at 25C	
B2	Blood agar base	Suspend 40g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Autoclave 15 min. at 121C Cool to 45 – 50C add 5 – 8% sterile Defibrinated Blood and Mix. Taking care to Avoid Bubble formation. Pour plates. Final PH 6.8 ± 0.2 at 25C	
B3	Brain Heart broth	Suspend 37g in 1L of Demin. Water. Dispense into smaller vessels. Autoclave 15 min. at 121C Final PH 7.4 ± 0.2 at 25C	
B4	BRILA broth	Suspend 40g in 1L of Demin. Water. Filled into tubes fitted with fermentation tubes. Autoclave 15 min. at 121C Final PH 7.2 ± 0.2 at 25C	BGB
C1	Cetrimid agar	Suspend 45.3g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Add 10 ml/L of Glycerol MERCK Cat No 1.04094 Autoclave 15 min. at 121C Final PH 7.2 ± 0.2 at 25C	

Code	Name	Preparation	Abb.
C2	Chapman agar	Suspend 146.5g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Autoclave 15 min. at 121C Final PH 7.0 ± 0.2 at 25C	
C3	Cooked meat broth	Place 1 tablet in a tube containing 10ml Demin. Water. Allow the tablet to disintegrate. Shake the suspension. Close the tubes with cotton wool plugs. Autoclave 15min. at 121C Final PH 7.4 ± 0.1 at 25C	
C4	Cooked meat medium	Suspend 12.5g in 100ml of Demin. Water. Mix thoroughly and allow to stand for 15 min. until all the particles are thoroughly wetted. Autoclave 15 min. at 121C & 15 lbs Final PH 7.2 ± 0.2 at 25C	
E1	Ec - Broth	Suspend 37g or 74g in 1L of Demin. Water. Dispense into tubes filled with fermentation tubes. Autoclave 15 min. at 121C Final PH 6.9 ± 0.2 at 25C	
E2	Endo agar	Suspend 39g in 1L of Demin. Water. by heating in a Boiling water bath or in a Current of Steam. Autoclave 15 min. at 121C A thick layer of medium. To deep a red color of medium. After solidification can be removed by adding a few drops (max 1ml/L) of a freshly prepared 10% solution of sodium sulfate and subsequent boiling. Final PH 7.4 ± 0.2 at 25C	

Code	Name	Preparation	Abb.
E3	Eosin methylene – blue Lactose Sucrose agar	Suspend 36g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Autoclave 15 min. at 121C Final PH 7.1 ± 0.2 at 25C	EMB
F1	Fluid lactose Medium	Suspend 13g in 1L of purified Demin. Water. Mix well and distribute into tubes with inverted Durham's tubes. Autoclave 15 min. at 121C & 15 lbs Final PH 6.9 ± 0.2 at 25C	
K1	Kliger iron agar	Suspend 55g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Dispense into tubes Autoclave 15 min. at 121C Final PH 7.4 ± 0.1 at 25C	
L1	Lactose broth	Suspend 13g in 1L of Demin. Water. Dispense into reagent tubes fitted with fermentation tubes Autoclave 15 min. at 121C Final PH 6.9 ± 0.2 at 25C	
L2	Lauryl sulfate agar	Suspend 92g in 1L of Demin. Water. Heat with frequent agitation until completely dissolved. Autoclave 15 min. at 121C Final PH 7.4 ± 0.2 at 25C	
L3	Lauryl sulfate broth	Suspend 36g in 1l of Demin. Water. Dispense into test tubes fitted with fermentation tubes. Autoclave 15 min. at 121C Final PH 6.8 ± 0.2 at 25C	LTB
L4	Litmus milk	Suspend 101g in 1L of Demin. Water Autoclave 15 min. at 121C & 15 lbs Final PH 6.8 ± 0.2 at 25C	

Code	Name	Preparation	Abb.
M1	Macconkey agar	Suspend 50g in 1L of Cold Distilled Water. Heat to boiling with frequent agitation. Autoclave 15 min. at 121C Cool to 45 – 50C. Mix well pour into sterile plates. Final PH 7.1 ± 0.2 at 25C	
M2	Methyl – red voges proskauer broth	Suspend 17g in 1L of Demin. Water. Dispense 5ml into tubes. Autoclave 15 min. at 121C Final PH 6.9 ± 0.2 at 25C	MR VP
N1	Nutrient agar	Suspend 20g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Autoclave 15 min. at 121C Final PH 7.0 ± 0.2 at 25C	NA
P1	Peptone water	Suspend 25.5g in 1L of Demin. Water. If desired dispense into smaller vessels. Autoclave 15 min. at 121C Final PH 7.2 ± 0.2 at 25C	
P2	Perfringens Selective agar	Suspend 40g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Autoclave 15 min. at 121C Final PH 7.0 ± 0.1 at 25C	SPS
P3	Plate count agar	Suspend 22.5g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Autoclave 15 min. at 121C Final PH 7.2 ± 0.2 at 25C	

R1	R- 2 A agar	Suspend 18.12g in 1L in Distilled Water. Heat to boiling and dissolve the medium completely. Autoclave 15 min. at 121C & 15 lbs (DO NOT OVERHEAT) Final PH 7.2 ± 0.2 at 25C	
----	-------------	---	--

Code	Name	Preparation	Abb.
S1	Sabouraud dextrose agar	Suspend 65g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Autoclave 15 min. at 121C (DO NOT OVERHEAT) Final PH 5.6 ± 0.2 at 25C	
S2	Selenite broth base	Dissolve 4g of sodium biselenite (L121) in 1L of Demin. Water then add 19g of CM 395 warm to dissolve. Mix well and fill out into containers to a depth of 15cm. sterilize in a boiling water bath or in a free flowing steam for 10 min. DO NOT AUTOCLAVE Final PH 7.1 ± 0.2 at 25C	
S3	SIM BIOS Medium	Suspend 30g in 1L of Cold Demin. Water. Heat to boiling with frequent agitation to dissolve completely. Distribute into into tubes. Autoclave 15 min. at 121C Final PH 7.3 ± 0.2 at 25C	
S4	Simmons citrate agar	Suspend 22.5g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Autoclave 15 min. at 121C Final PH 6.6 ± 0.2 at 25C	
S5	SS agar	Completely Suspend 60g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. DO NOT AUTOCLAVE Cool Rabitly and Pour Plates to give a thick layer of Medium. Final PH 7.0 ± 0.2 at 25C	

Code	Name	Preparation	Abb.
T1	TCBS	Completely Suspend 88g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. DO NOT AUTOCLAVE Final PH 8.6 ± 0.2 at 25C	
T2	Triple sugar iron agar	Suspend 65g in 1L of Demin. Water by heating in a Boiling water bath or in a Current of Steam. Dispense into tubes. Autoclave 15 min. at 121C Allow to solidfy to give agar slants. Final PH 7.4 ± 0.2 at 25C	TSI
Y1	Yeast extract granulated	-	
Z1	Zhelaten powder	-	